

Homogenizer의 단짝친구, **Bead**

우리 **Bead**는요,
크기도 재질도 다양한 **Bead**,
고르기 고민되실 때는 바이오코드에
상담을 요청하시면,
샘플에 따라 추천해드립니다!

Bead	
BCS-BT13721	Pre-filled tube kits I, 80 prep (Lung, Liver 등 Soft tissue 에 최적화) (3 mm zirconium bead 7ea + 5 mm stainless bead 1ea/tube)
BCS-BT13722	Pre-filled tube kits II, 80 prep (Heart, tail 등 Hard tissue 에 최적화) (3 mm zirconium bead 7ea + 5 mm stainless bead 2ea/tube)
BCS-BT30302	Stainless bead (3 mm), 200 g
BCS-BT30304	Stainless bead (3 mm), 400 g
BCS-BT30505	Stainless bead (5 mm), 500 g
BCS-BT31302	Stainless bead (3 mm), sterile, 200 g
BCS-BT31304	Stainless bead (3 mm), sterile, 400 g
BCS-BT31505	Stainless bead (5 mm), sterile, 500 g
BCS-BT40011	Zirconium bead (0.1 mm), 100 g
BCS-BT40055	Zirconium bead (0.5 mm), 500 g
BCS-BT40105	Zirconium bead (1 mm), 500 g
BCS-BT40305	Zirconium bead (3 mm), 500 g
BCS-BT90200	2 mL High-qual, Screw tube and cap, 500/pk

Bead Selection Guide

		Zirconium				Stainless steel
		0.1 mm	0.5 mm	1 mm	3 mm	3 mm
Tissue	Skin				●	
	Hair				●	●
	Nail				●	●
	Bone				●	●
	Teeth				●	●
	Lung				●	
	Muscle				●	●
	Liver				●	
	Brain				●	
	Heart				●	
	Kidney				●	
	Pancreas				●	
	Fly				●	●
	Tick				●	●
	Gonad				●	
	Artery				●	
	Intestine				●	
	Cornea				●	●
	Tail				●	●
	Spleen				●	
	Ear				●	
Microbial	Gram +	●	●			
	Gram -	●	●			
	Yeast		●	●		
	Fungi		●	●		
	Bacteria	●	●			
	Spores		●	●		
	Mold		●	●		
Plant	Soft plant tissue					●
	Tough plant tissue					●
	Leaves					●
	Arabidopsis					●
	Seeds					●
	Corn					●
	Nuts					●
	Rice					●
	Wheat					●
	Stems					●

Plant의 경우, 3 mm Stainless steel bead와 5 mm Stainless steel bead를 혼합하여 사용하면 파쇄 효율을 더욱 높일 수 있습니다.